## PI800 ILLUMINATOR

ILLUMINATEUR PL800 • PL800 ILLUMINATOREN • ILUMINADOR PL800

A PRODUCT OF DOLAN-JENNER INDUSTRIES

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### OPERATION MANUAL

#### Introduction

The Fiber-Lite PL800 Series is a family of 150 Watt quartz halogen, fiber optic illuminators.

#### INITIAL SETUP

- Remove the illuminator from the carton.
- Loosen the fiber optic nosepiece thumbscrew.
- Insert the fiber optic into the illuminator nosepiece.
- Tighten the thumbscrew by hand to make a secure connection to the fiber optic. *Use of pliers or other tools is not recommended.*
- Insert the AC power cord into the power entry at the rear of the illuminator. *Use only approved power cord supplied with the illuminator.*

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## FIBER OPTIC CONNECTION

Two nosepiece styles are available with the PL800 series of fiber optic illuminators. A-type with a 0.985 in. (25mm) ID, or B-type with a 0.590 in. (15mm) ID smaller than the 0.985 in. ID.

A-type nosepieces require the use of adapters to mate fiber optic light guides smaller than the 0.985 in. ID. Adapters fit into the nosepiece and reduce the ID to accept the chosen fiber optic. The fiber optic and the adapter are secured with the thumbscrew. The following A-type adapters (SX-5, SX-6 & SX-7) are made with an idler which allows quick disconnect of the fiber optic.

B-Type nosepieces will accept self-supporting goosenecks (EEG2823C, EEG3922C, BG2820C, etc.) and ringlight (A3739B) fiber optics without the need for an adapter. Fiber optics smaller than the 0.590 in. (15mm) ID will require the use of B-type adapters (SX-5B, SX-6B & SX-7B).

## OPERATION

All the PL800 series illuminators are equipped with a front panel manual solid state intensity control. Some models also have an external interface (9-pin, D style connector) for remote intensity control via a 0-5 VDC signal on the back panel.

First insert the power cord into the power entry module at the rear of the illuminator. Next, insert the power cord into a 3 wire grounded AC power receptacle.

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**MANUAL:** Press the ON(1)/OFF(0) rocker switch on the front panel to ON(1). The illuminator will light and is ready for service. The illuminator intensity is controlled by the rotary control located on the front panel of the illuminator. The 0 position (when control is turned fully counter-clockwise) corresponds to the lowest illuminator intensity. The 100 position (when control is turned fully clockwise) corresponds to the highest illuminator intensity. *NOTE: Continuous operation of the illuminator at the highest intensity level will yield rated lamp life. Operating the illuminator at reduced intensity can result in significantly extended lamp life.*

**REMOTE:** Switch Manual/Remote Switch on the back panel to the Remote position. Intensity will now only be controlled via the 0-5 VDC analog input signal. Press the front panel ON(1)/OFF(0) rocker switch to ON(1). The illuminator will light and is ready for service. *NOTE: The front panel intensity control is deactivated in remote mode.*

The 0-5 VDC analog input signal interface is the 9-Pin connector located at the rear of the illuminator. See diagram for Pinout.

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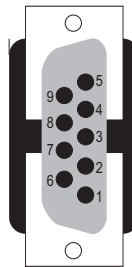
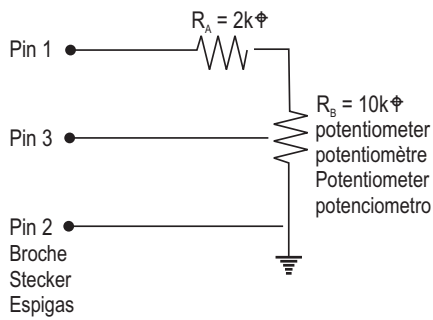
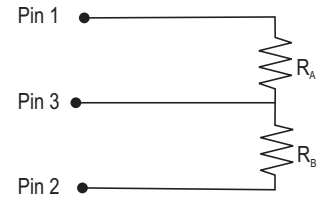
II

I. The +6 VDC at 5mA max. signal available at Pin #1 can be used as a voltage source for a remote potentiometer control (use  $R_A \geq 2k\ \Omega$  minimum resistance).

II. For fixed intensity use  $\geq 2k\ \Omega$  resistance. ( $R_A + R_B \geq 2k\ \Omega$ )

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**⚠ WARNING • AVERTISSEMENT • WARNUNG • PRECAUCIÓN ⚠**

Risk of electrical shock. Remove power plug before lamp replacement and wait for hot lamp to cool.

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**LAMP MODULE REPLACEMENT**

1. Turn the illuminator intensity control fully counterclockwise (the 0 position) and run the illuminator with the fan for several minutes. Wait until the nosepiece is cool to the touch. Press the ON(1)/OFF(0) rocker switch to the OFF(0) position.
2. Remove the AC line cord from the AC power receptacle.
3. Release lamp module from the power supply by turning both retaining screws counterclockwise using a straight blade screwdriver. *NOTE: Both retaining screws will be disengaged from the power supply, but will remain in place in the lamp module.*
4. Remove the lamp module from the power supply by slowly pulling outward.

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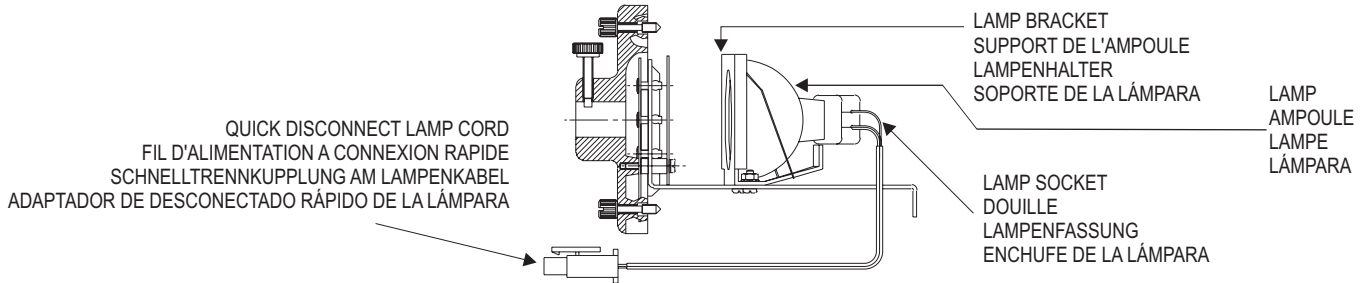
5. Release the lamp module by unlatching the quick disconnect lamp cord. See diagram III.
6. Attach new lamp module by connecting the quick disconnect lamp cord.
7. Insert lamp module into power supply. Make sure lamp cord does not interfere.
8. Secure lamp module by tightening the (2) retaining screws clockwise.
9. Reattach AC line cord and the illuminator is ready for service.

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**III LAMP MODULE • MODULE DE LA LAMPE • LAMPENMODULS • MÓDULO DE LA LÁMPARA**

**LAMP REPLACEMENT**

1. Follow steps 1-4 of Lamp Module Replacement.
2. Check the lamp assembly to verify that the lamp and socket are cool before proceeding.
3. Lift and remove the lamp from the lamp holder by grasping the rear of the lamp adjacent to the lamp socket.
4. Remove the lamp from the socket by holding the lamp socket and gently pulling the lamp reflector.
5. Discard the lamp.
6. Insert the replacement lamp into the lamp socket. *CAUTION: Do not touch the interior of the lamp reflector, the lamp envelope or the lamp pins with your fingers. Touching the interior of the lamp reflector, the lamp envelope or the lamp pins will result in significant shortening of the lamp life. Handle the lamp only by the exterior of the reflector or the area adjacent to the pins.*

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**WARNING**

Replace the fuse with the correctly rated fuse as listed on the label on the back of the illuminator. Use of an improper fuse can create a hazardous situation.

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7. Insert the replacement lamp and socket into the lamp holder.
8. Follow steps 7-9 of **Lamp Module Replacement**.

**FUSE REPLACEMENT**

1. Press the ON(1)/OFF(0) switch to the OFF(0) position.
2. Remove the AC line cord from the AC power receptacle.
3. Remove the AC line cord from the power entry module at the rear of the illuminator. The fuse drawer is part of the power entry module. The drawer is located directly beneath where the AC line cord plugs in.
4. **For 115 VAC illuminators:** Pull out the fuse drawer. Remove the blown fuse that is positioned closest to the illuminator and discard. The second fuse is the spare.
5. Place the replacement fuse into the fuse drawer. The fuse will work in either orientation.
6. Push the fuse drawer until it "clicks" into position.
7. **For 230 VAC illuminators:** Push down the center latch until it locks in the down position. Remove the fuse drawer. Two fuses are required.
8. Remove the blown fuse(s) from the drawer and discard.

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9. Place the replacement fuse(s) into the fuse holder by inserting either end of the fuse into the fuse holder.

10. Push the fuse drawer until it "clicks" into position.

11. Attach the AC line cord to power entry module at the rear of the illuminator. The illuminator is now ready for service.

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*Only above parts are replaceable. Return illuminator to factory for warranty service. Attempts to replace other parts voids warranty.*

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## PERFORMANCE STATEMENT

Dolan-Jenner Industries, Inc. (DJI) recognizes that its illuminator products may be used under an almost unlimited variety of conditions. As such, we are prepared to assist the customer in the selection and application of any of these products. This includes application engineering, sample testing and other means as determined by DJI.

Where DJI has made specific recommendations for its products, systems, or detection techniques (based on complete and detailed information furnished by the customer) we will extend every effort to assure that the customer is satisfied with the performance of our products. Continual development and improvement of DJI products may require changes in details that do not coincide with descriptions or illustrations shown. All fiber optic bundle diameters are nominal.

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## WARRANTY ON LIGHT SOURCES

Dolan-Jenner Industries, Inc. (DJI) warrants its products to be free from defective material and workmanship. Any light source or parts thereof which are determined by DJI to be defective within two years from shipment date will be replaced or repaired at our option. This policy is effective Aug. 1, 2005 and is not retroactive. All fiber optics are warranted for one year. This warranty specifically excludes both incandescent and quartz-halogen lamps, and optical filters.

Any products which, in our opinion, have been subjected to misuse, incorrect wiring, or where installation procedures are not in accordance with the instruction manual, are excluded from this warranty. Nor does this warranty extend to products on which repairs or alterations have been made outside the factory, or on which the identification or serial number has been altered or to accessories not of our manufacture.

Our obligation with respect to products or parts covered by this warranty shall be limited to repair or replacement, F.O.B., Boxborough, Massachusetts. In no event shall DJI be held liable for consequential or special damages, or for transportation, installation, adjustment, or other expenses which may arise in connection with such products or parts. This warranty is in lieu of all other statements or warranties or guarantees, written or implied, by DJI or its authorized representatives.

**Important: Please contact factory for a return authorization number prior to shipping merchandise to factory.**

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**TECHNICAL DATA**

Lamp	150 W Quartz halogen, 21V (EKE, EJV or EJA0 or 20V (DDL) as ordered
Voltages	115 VAC, 200 W, 50/60 Hz or 230 VAC, 1 A, 50/60 Hz (Set at factory)
Lamp Life	200-10,000 Hours depending on intensity
Fuse	115 VAC: 3 Amp, 250V, time delay, 5 x 20 mm 230 VAC: 3 Amp, 250V, time delay, 5 x 20 mm
Color Temperature	3100-3400 degrees Kelvin depending on lamp used
Noise Level	Approximately 21 dB (A)
Safety Approvals	UL
Dimensions	H x W x D (inch): 4.6 x 7.25 x 9.5 H x W x D (cm): 11.7 x 18.4 x 24.3
Weight	12 lbs. (5.5 kg.)
Max. Housing Temp.	9°C above ambient

**Environmental Conditions Use Specifications**

Type of Use	Indoor at max. altitude of 2,000 M
Installation Category	II
Temperature Range	5°C to 40°C
Relative Humidity	@ 80% maximum up to 31° C decreasing linearly to 50% relative humidity @ 40° C
Pollution Degree	2

**NOTICE: if the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.**

**Corporate Headquarters and Factory:**

Dolan-Jenner Industries, Inc.  
159 Swanson Rd.Boxborough, MA 01719, USA  
Toll Free: 1.800.83.FIBER  
Telephone: 978.263.1400  
Fax: 978. 264.0292

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